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1 Safety notes

Protect the sensitive bronze band of the torsion pendulum from uncontrolled mechanical loading:

- Do not unscrew the oscillating system's locking screws (7) - see Fig. 2 - before the device has been correctly assembled and brought into position.
- Always lock the oscillating system when the device is not in use. In particular, make sure it is locked during transport and assembly.

The knurled screw (3.1) - see Fig. 2 - for fixing the torsion head (preadjusted on delivery) is only to be slightly loosened if a fine adjustment of the zero point proves necessary on putting the correctly-assembled torsion balance into operation. The grub screw (3.2) - see Fig. 2 - for fixing the pendulum holder is only to be loosened when replacing the torsion band as described in section 4.

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2 Description, technical data, scope of supply (see Fig. 2)

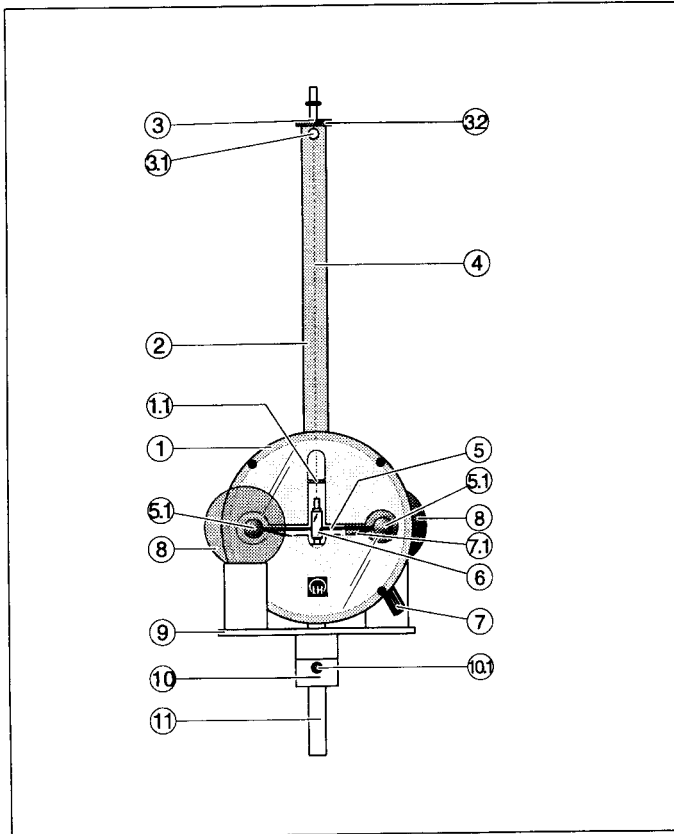
Fig.2
Gravitations-Drehwaage.
Gravitation torsion balance

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- ① Metal housing (15 cm dia.), with metal ring and glass covers; two-part slide (1.1) to prevent interference from convection currents between the housing and protective tube (2)
- ② Protective tube (25 cm long) for torsion band (4)
- ③ Torsion head with pendulum holder, can be rotated for adjustment of the pendulum equilibrium position when screw has been loosened (3.1); pendulum holder fixed with grub screw (3.2)
Important!
Do not loosen the knurled screw (3.1) unless it is essential to carry out a zero-point adjustment in accordance with section 3.5. Only loosen the grub screw (3.2) when the torsion band is being replaced as described in section 4.
- ④ Torsion band made from bronze, 26 cm long
Spare part no.: 683 21
- ⑤ Dumbbell-shaped pendulum body, consisting of 2 lead balls (5.1) on a metal rod