



- \* Microphone: low impedance, 50-15,000 Hz frequency response (replacement battery = 1.5 V Eveready A76 or equivalent)

### Additional Equipment Needed:

- \*3. PASCO Microwave Transmitter
- \*4. PASCO Microwave Receiver
- \*5. 9 V Power Adapter
- 6. an oscilloscope, or an audio amplifier and speaker, to monitor the output.
- 7. Banana plug connectors. (between the oscilloscope and the modulation)

\* Items 3-5 are provided in the PASCO Microwave Optics System (WA-9314B or WA-9316)

### Operation

Set up the equipment as shown in the preceding illustration. Turn on the microphone with the ON/OFF switch found on the microphone battery compartment. Point the Microwave Transmitter toward the Microwave Receiver. Then:

- To monitor the modulating signal, plug your oscilloscope into the banana plug connectors labeled SCOPE on the Microwave Modulator. (see Figure)
- To monitor the signal at the receiving end, plug your oscilloscope or audio amplifier w/speaker (not included) into the banana plug connectors labeled OUTPUT on the top of the Microwave Receiver. (see Figure)
- If you wish to modulate the microwaves with the triangular wave signal, unplug the microphone. Adjust the frequency, amplitude, and bias of the modulating signal using the controls on the Microwave Modulator. (The BIAS control adjusts the DC voltage level of the Gunn Diode in the Microwave Transmitter. The modulating signal then varies the Gunn Diode voltage about this DC voltage. Proper adjustment of the BIAS is necessary to produce a linear, unclipped microwave signal from the Gunn diode.)

## Experiments

### 1. ADDING AN AUDIBLE OUTPUT TO MICROWAVE EXPERIMENTS


The impact of most microwave experiments can be improved by adding an audible output. The maxima and minima of an interference pattern, for example, seem somehow more real when they can be heard in the variations in loudness of an audible tone, instead of just seen as fluctuations in a meter reading.

Any of the experiments in the manual that accompanied your PASCO Microwave Optics System can be performed with an audible output. Just follow the instructions listed above for using the triangular wave output of the Modulator with an audio amplifier and speaker.


### 2. MICROWAVE COMMUNICATIONS

With the simultaneous use of an oscilloscope and an audio amplifier and speaker, this system is excellent for demonstrating microwave communications. The input signal can be the triangular wave output of the modulator, or it can come from the microphone or even a tape player.

#### Typical View of Dual Trace Oscilloscope

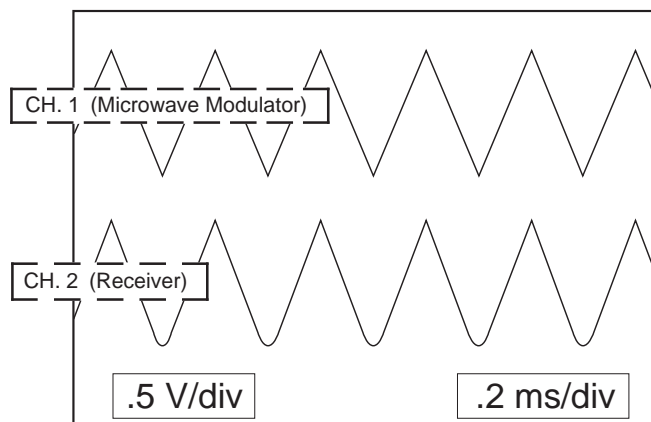


#### Typical Equipment Settings

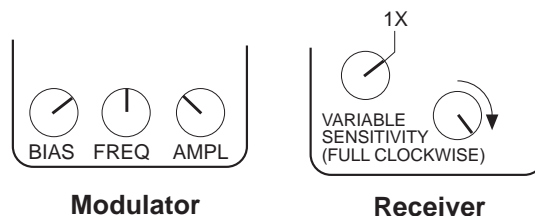

Modulator

Receiver

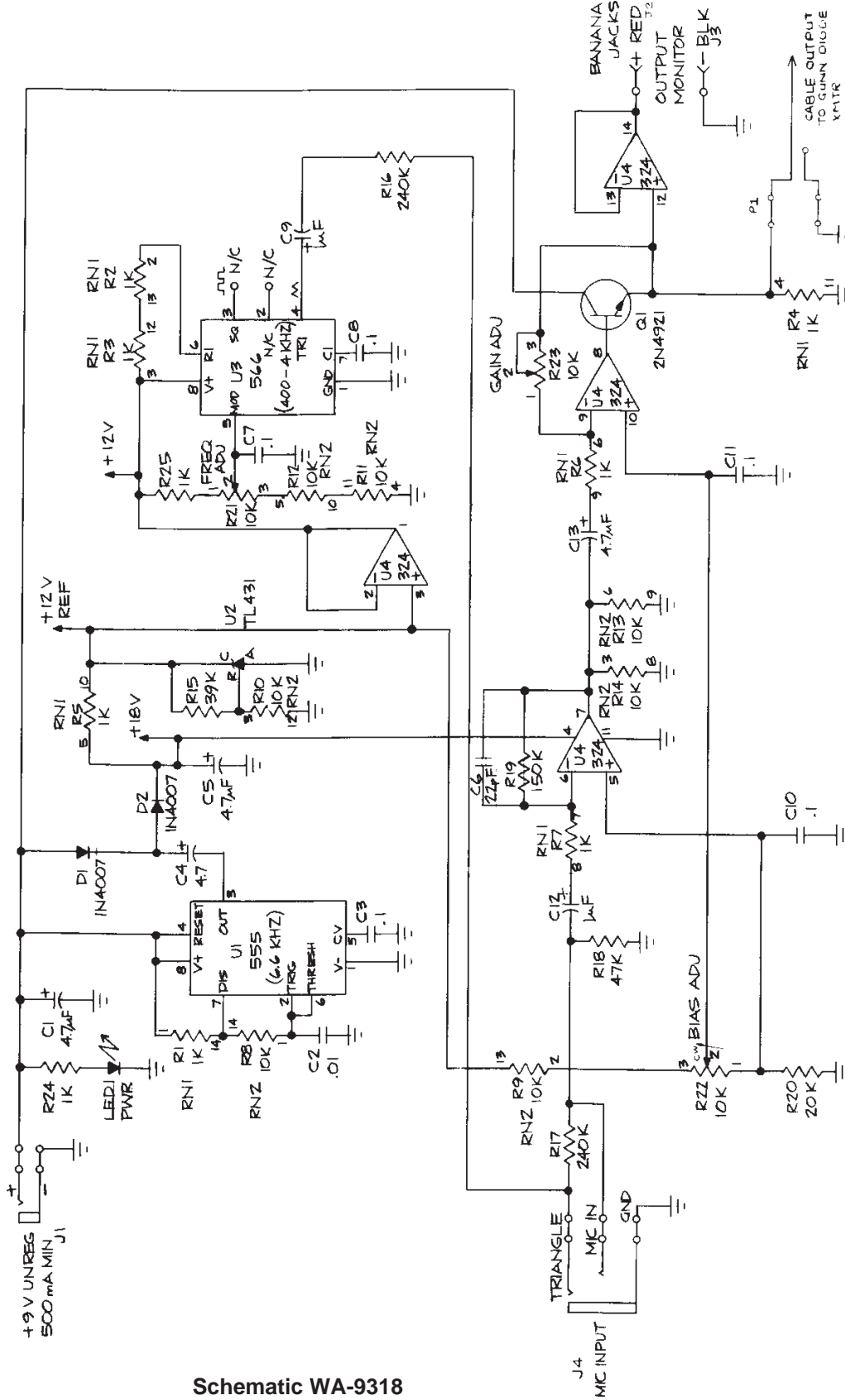

Schematic WA-9318  
(Drawing No.956-02857)

## Limited Warranty

PASCO scientific warrants this product to be free from defects in materials and workmanship for a period of one year from the date of shipment to the customer. PASCO will repair or replace, at its option, any part of the product which is deemed to be defective in material or workmanship. This warranty does not cover damage to the product caused by abuse or improper use. Determination of whether a product failure is the result of a manufacturing defect or improper use by the customer shall be made solely by PASCO scientific. Responsibility for the return of equipment for warranty repair belongs to the customer. Equipment must be properly packed to prevent damage and shipped postage or freight prepaid. (Damage caused by improper packing of the equipment for return shipment will not be covered by the warranty.) Shipping costs for returning the equipment, after repair, will be paid by PASCO scientific.

## Equipment Return

Should the product have to be returned to PASCO scientific for any reason, notify PASCO scientific by letter, phone, or fax BEFORE returning the product. Upon notification, the return authorization and shipping instructions will be promptly issued.

► **NOTE: NO EQUIPMENT WILL BE ACCEPTED FOR RETURN WITHOUT AN AUTHORIZATION FROM PASCO.**

When returning equipment for repair, the units must be packed properly. Carriers will not accept responsibility for damage caused by improper packing. To be certain the unit will not be damaged in shipment, observe the following rules:

- ① The packing carton must be strong enough for the item shipped.
- ② Make certain there are at least two inches of packing material between any point on the apparatus and the inside walls of the carton.
- ③ Make certain that the packing material cannot shift in the box or become compressed, allowing the instrument come in contact with the packing carton.

Address: PASCO scientific  
10101 Foothills Blvd.  
Roseville, CA 95747-7100

Phone: (916) 786-3800  
FAX: (916) 786-3292  
email: techsupp@pasco.com  
web: www.pasco.com